Levenhuk DTX

Digital

EN	User I
CZ	Návod
DE	Bedie
ES	Guía d
PL	Instru
RU	Инстр

500x ZOOM	5
ION BATTERY	m

leve

Levenhuk DTX 500 LCD


Digital Microscope

EN

Never, under any circumstances, look directly at the light source or another bright source of light or at the sun, as this may cause PERMANENT damage to your eyes and may lead to BLINDNESS.

Microscope parts

- | | |
|---------------------------|-----------------------------------|
| 1 LCD display | 9 Brightness adjustment wheel |
| 2 MicroSD card slot | 10 Magnification adjustment wheel |
| 3 Fine focus knob | 11 Power cable socket |
| 4 Microscope optical tube | 12 RCA socket |
| 5 Stage | 13 USB socket |
| 6 Coarse focus knob | 14 Battery compartment |
| 7 Focus lock knob | 15 Battery status indicator |
| 8 Shutter button | 16 Function keys |

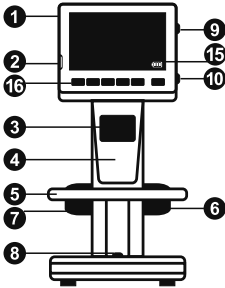The kit includes: microscope, rechargeable battery, power cable, USB cable, RCA cable, Levenhuk DTX installation CD, cleaning cloth, user guide, calibration scale.

Microscope assembly

Open the battery compartment at the bottom of the microscope base. Insert the battery into the battery compartment matching the indicated polarity. The battery has to be fully charged before turning the microscope on for the first time. Plug the power cable into the power cable socket on the microscope body (11) and connect it to a power source. The battery status indicator (15) will turn red. Once the battery is fully charged, the indicator will go out. You can now disconnect the microscope from the power source and turn it on. The fully-charged battery allows you to use the microscope for two hours. The illumination will grow dim and the image quality will decrease to inform you that the battery is low. Charge the battery, as described above.

A MicroSD card (purchased separately) is required to save captured images and recorded videos. Turn the microscope off and insert the MicroSD card into the corresponding slot on the microscope body until it locks in place with a click. Do not apply excessive force when inserting the MicroSD card. Try rotating the card, if it does not lock in place with ease. Format the card before using it.

To turn the microscope on press the power button (16). To turn the microscope off press the power button (16) again (refer to Power saving for more information).

Image cap

Sequence

00013/01618

Captured
of images
(based on

MicroSD s

5.0M

Current im

Battery st

Note: inte
the magn

Using the microscope

Capturing an image

Turn the microscope on and place a specimen on the stage. Adjust the illumination brightness, so that the specimen is evenly lit. Change magnification and focus your view, if necessary, then lock the results in place with the focus lock knob (7). Press the shutter button (8).

Recording a video

Press the ◀ key (16) to switch to video mode. The icon in the upper left corner of the screen will change in confirmation. Press the shutter button (8) to start recording and press it again to stop.

Playing recorded videos

Press the ◀ key (16) to switch to playback mode. The icon in the upper left corner of the screen will change in confirmation. Select the desired video clip with scroll keys (16) and press **OK** to play it.

Deleting files

When browsing saved files, press the **Wrench** key (16), select **Delete...** press **OK** to confirm selection (you can choose to delete everything or just the selected file from the displayed menu).

Protecting files

When browsing saved files, press the **Wrench** key (16), select **Protect** press **OK** confirm selection (you can choose to protect everything or just the selected file from the displayed menu).

Slideshow

When browsing saved files, press the **Wrench** key (16), select **Slide Show** and press **OK** to confirm selection.

Interface color mode

When browsing saved files, press the ▲ key to switch between existing interface color modes.

Magnification

Scroll the magnification adjustment wheel (10) to zoom your view in or out.

Note: magnification range of Levenhuk DTX 500 LCD is 1x–4x.

Transferring files to a computer

Connect the microscope to your computer with a USB cable. When the connection appears, choose to use the microscope as an external drive.

Main menu

Press the **Wrench** key (16) to open the main menu. Select **Resolution** to adjust resolution, adjust image quality and sharpness settings, and press **OK** to capture.

Resolution

Select **Resolution** from the main menu choose the desired resolution and press **OK** to confirm selection.

Note: model supports the following resolutions

Sequence capture

Select **Timer** from the main menu and set the desired interval. Press **OK** to open the **Frames** dialog window. Confirm your selection. The **Frames** dialog will appear where you can set the interval between captures. Press **OK** to confirm and ◀ to return to capture mode. Press **OK** to start capture. The icon in the upper left corner will change in confirmation. When a frame has been captured, the screen will return to stop sequence capture. Press **OK** at any time to stop sequence capture.

Quality

Select **Quality** from the main menu and choose the desired quality. Press **OK** to confirm the compression rate of captured images.

Sharpness

Select **Sharpness** from the main menu and choose the desired sharpness. Press **OK** to confirm selection.

Language

Select **Language** from the main menu and choose the desired language. Press **OK** to confirm selection.

Power saving

Select **Power Save** from the main menu and choose the desired power saving mode. Press **OK** to confirm. The microscope automatically shuts down.

Display frequency

Select **Frequency** from the main menu and choose the desired display frequency. Press **OK** to confirm selection.

Date and time display

Select **Date Stamp** from the main menu and choose one of the existing settings. You can opt to display the date or both date and time. Press **OK** to confirm selection.

TV connection

Select **TV Out** from the main menu and choose one of the existing settings. Press **OK** to confirm selection.

Date and time

Select **Date/Time** from the main menu and set the current date and time. Press **OK** to confirm entry. You can also choose one of the following date formats: YY/MM/DD, DD/MM/YY or MM/DD/YY.

Format SD card

Select **Format SD** from the main menu to format the installed memory card.

Settings reset

Select **Sys. Reset** from the main menu to revert back to factory settings.

Hardware version

Select **Version** from the main menu to check the current hardware version.

Connecting the microscope to a TV-set

Plug the RCA cable into the microscope and connect it to your TV-set (refer to the User Guide of your TV-set for additional information). Turn on the TV and select the correct channel from the corresponding menu of the TV-set (refer to the User Guide of your TV-set for additional information).

PortableCapture Pro software

Insert the installation CD into your CD drive. The Main Menu window will be displayed. Select **PortableCapture software** to install the PortableCapture Pro application. Keep in mind that a system restart will be required during installation. The installation CD contains a copy of this User Guide, which you can read by clicking the corresponding entry in the Main Menu.

Connect the microscope to your PC with a USB cable and run the PortableCapture Pro application. A pop-up message will warn you if the microscope is not connected to the PC.

Note: function keys are disabled when the microscope is connected to a computer as a camera. The microscope is operated through the PortableCapture Pro application. All the images and videos will be saved to the computer's hard drive.

Capture an image

Start and stop video recording

Sequence capture (you can capture images or videos, interval between images or videos to be captured)

Full screen mode. To exit full screen mode, press F11 on your keyboard or double-click the mouse button

Save as

Copy to clipboard

Copy file

Previous image

Next image

Undo the last action

Repeat the last action

File

Photos Directory: set a directory for captured images.

Videos Directory: set a directory for recorded videos.

Options

Resolution: set image resolution.

Date/Time: hide or display date and time of capture during preview.

Language: change the language of the user interface.

Crossing: hide or display a grid on images.

Full screen mode: go full screen.

Capture

Photo: capture an image.

Video: record a video.

Capturing an image

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu.

Recording a video

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu.

Sequence capture

Click the corresponding icon on the toolbar. A dialog window will appear where you can set a number of parameters for sequence capture. Click the **Photo** radio button to set the time to begin capturing images, interval between them and the total number of images. Click the **Video** radio button to set the time to begin recording videos, their duration, interval between them and the total number of clips.

Calibrating

Before calibrating the microscope, turn the grid display on. To do so, select **on** from **Options > Crossing**. By default, the grid is visible. Place the calibration scale on the stage and focus your view. Make sure that the vertical axis of the grid is parallel to scale divisions and capture an image. Double-click the image thumbnail to open it in a separate window.

Click the corresponding icon on the toolbar to begin calibration. Using your mouse, choose two points on the image (you have to know the actual distance between these two points). Once you choose the second point, a dialog window will open, where you have to enter the known distance in the **Actual dimension** field. The application will automatically calculate the magnification of the image. Click **OK** to close the dialog window. The calculated magnification will be displayed in the **Magnification** field.

Note: chosen points should form a horizontal line.

You can check calibration results by measuring with the application's ruler tool. Click the measurements icon, choose the tool, place it on the image. If the measured distance is equal to the known distance, the calibration is successful.

Recalibration

The calibration process has to be repeated if the magnification changes during observations.

Click the corresponding icon on the toolbar to begin recalibration. Choose the tool from **Reset magnification**. Repeat the calibration process: place the tool on the image, magnification and capture a few images. Double-click the image thumbnail to open the resulting image in a preview window. Click the corresponding icon on the toolbar to **picture magnification** from **Set magnification**. A dialog window that appears and click **OK**.

If you want to use the same magnification when recalibrating, click the corresponding calibration icon on the toolbar and select **Set magnification**.

Measurements

It is recommended to calibrate the system before measurements.

Any Angle Line. Press and hold the left mouse button, move the mouse to draw a line. Release the mouse button to complete the line.

Continuous Line. Use this tool to draw a curve.

Radius Circle. Draw a straight line on the image. Release the mouse button to draw the radius, as well as circumference and area of the circle.

Diameter Circle. Draw a straight line on the image. Release the mouse button to calculate the diameter, as well as circumference and area of the circle.

Three Points Angle. Place three points on the image. Release the mouse button to automatically calculate the angle value.

Draw and text

PortableCapture Pro allows you to add drawing tools to the image. Click the corresponding icon on the toolbar to draw a line, rectangle, circle, etc. Click the **Pen** icon on the toolbar to draw text. Click the **Pen** icon from the drop-down menu and draw something on the image. In the preview window, you can edit font type and color to be used.

Specifications

Image sensor	5Mpx (up to 12Mpx by interpolation)
Magnification	20x–200x–500x
Available image resolutions	12Mpx, 8Mpx, 5Mpx, 3Mpx, 1.3Mpx, VGA
Photo/video format	*.jpeg/*.avi
Memory card	microSD up to 32Gb (not included)
Objective lens	High-quality lens
Display	3.5" TFT-display, 4:3 ratio
Illumination	8-LED system with variable brightness
Power source	Li-ion battery (3.7V/1050 mAh)
	Operation time: 2 hours, charging time: 2 hours
Frame rate	30fps
Power cable	100–240V, 50/60Hz; 5V, 1A output
OS language	English, German, French, Spanish, Italian, Portuguese, Japanese, Chinese
Software language	English, German, French, Spanish, Russian, Italian, Portuguese, Polish, Dutch, Japanese, Korean, Chinese
Dimensions (HxWxD)	227 x 104 x 149mm
Weight	760g (with the battery)

Levenhuk reserves the right to modify or discontinue any product without prior notice.

System requirements

Operating system: Windows XP SP2/Vista/7/8, Mac 10.6–10.8,
CPU: at least P4 1.8GHz, RAM: 512MB, interface: USB 2.0

Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a lifetime warranty against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk. This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

For further details, please visit our web site: www.levenhuk.com/warranty

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Levenhuk® is a registered trademark of Levenhuk, Inc. © 2006-2015 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 924-D East 124th Ave. Tampa, FL 33612 USA