# Levenhuk DTX

## Digital

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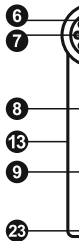
## Digital Microscope

EN

Never, under any circumstances, look directly at the sun or another bright source of light or use the microscope as a telescope, as this may cause PERMANENT damage to your eyes and may lead to BLINDNESS.

### Microscope parts

- |                         |  |                             |
|-------------------------|--|-----------------------------|
| 1 Shutter button        | 9 MicroSD card slot                    | 17 OK                       |
| 2 Zoom adjustment wheel | 10 Wrist strap clip                    | 18 ▼ (DOWN)                 |
| 3 RCA socket            | 11 Focus wheel                         | 19 ◀ (LEFT/ BACK)           |
| 4 USB socket            | 12 Brightness adjustment wheel         | 20 Power button             |
| 5 Power cable socket    | 13 Battery compartment                 | 21 Speaker                  |
| 6 LEDs                  | 14 LCD display                         | 22 Microphone               |
| 7 Objective lens        | 15 ▲ (color mode switch and scrolling) | 23 Battery status indicator |
| 8 Tripod threaded hole  | 16 Menu                                |                             |

The kit includes: microscope, tripod, rechargeable battery, power cable, USB cable, RCA cable, Levenhuk DTX installation CD, pouch, wrist strap, cleaning cloth, user guide.

### Microscope assembly

Open the battery compartment and insert the battery into the battery compartment matching the indicated polarity. The battery has to be fully charged before turning the microscope on for the first time. Plug the power cable into the power cable socket on the microscope body (5) and connect it to a power source. The battery status indicator (23) will turn red. Once the battery is fully charged, the indicator will go out. You can now disconnect the microscope from the power source and turn it on. The fully-charged battery allows you to use the microscope for two hours. The illumination will grow dim and the image quality will decrease to inform you that the battery is low. Charge the battery, as described above.

A MicroSD card (purchased separately) is required to save captured images and recorded videos. Turn the microscope off and insert the MicroSD card into the corresponding slot on the microscope body until it locks in place with a click. Do not apply excessive force when inserting the MicroSD card. Try rotating the card, if it does not lock in place with ease. Format the card before using it.

To turn the microscope on, open the LCD display or press the power button (20). To turn the microscope off, fold the LCD display or press the power button (20) again (refer to Power saving section for more information).

Image cap

Sequence

00013/01618

Captured  
of images  
(based on

5.0M

MicroSD s

Current im

Battery st

Note: inte  
the magn

## Using the microscope

## Capturing an image

Turn the microscope on and point it at an object. Make sure the LCD display is fully unfolded. Adjust the illumination brightness, so that the specimen is evenly lit. Change magnification, then focus your view and press the shutter button (1) to capture an image.

## Recording a video

Press the ◀ key (19) to switch to video mode. The icon in the upper left corner of the screen will change in confirmation. Press the shutter button (1) to start recording and press it again to stop.

### Playing recorded videos

Press the **◀** key (19) to switch to playback mode. The icon in the upper left corner of the screen will change in confirmation. Select the desired video clip with scroll keys and press **OK** to play it.

## Deleting files

When browsing saved files, press the **Wrench** key (16), select **Delete...** press **OK** to confirm selection (you can choose to delete everything or just the selected file from the displayed menu).

## Protecting files

When browsing saved files, press the **Wrench** key (16), select **Protect** press **OK** confirm selection (you can choose to protect everything or just the selected file from the displayed menu).

## Slideshow

When browsing saved files, press the **Wrench** key (16), select **Slide Show** and press **OK** to confirm selection.

### Interface color mode

When browsing saved files, press the ▲ (15) key to switch between existing interface color modes.

## Zoom

Scroll the zoom adjustment wheel (2) to zoom your view in or out.

**Note:** zoom range of Levenhuk DTX 500 Mobi is 1x–4x.

## Attaching the tripod

Place the tripod on a stable surface. Connect the tripod to the threaded hole (10) on the microscope and rotate the microscope until it locks in place. Do not apply excessive pressure during rotation, as this may damage both the microscope and the tripod. Tripod height may be adjusted, if necessary.

## Transferring files to a computer

Connect the microscope to your computer with the cable that appears. choose to use the microscope as an external device.

**Main menu**

Press the **Wrench** key (16) to open the main menu for resolution, adjust image quality and sharpness, and capture.

## Resolution

Select **Resolution** the main menu choose the Press **OK** to confirm selection.

*Note: model supports the following resolutions*

### Sequence capture

**Select Timer** from the main menu and set the **Frames** dialog window. Confirm your selection appear where you can set the interval between confirm and ◀ to return to capture mode. Press The icon in the upper left corner will change in has been captured, the screen will return to st at any time to stop sequence capture.

## Quality

Select **Quality** from the main menu and choose the compression rate of captured images.

### Sharpness

Select **Sharpness** from the main menu and ch

## Language

Select **Language** from the main menu and choose the language you want to use to confirm selection.

### Power saving

Select **Power Save** from the main menu and select **On**. The system automatically shuts down.

### Display frequency

Select **Frequency** from the main menu and choose **OK** to confirm selection.

### Date and time display

Select **Date Stamp** from the main menu and choose one of the existing settings. You can opt to display the date or both date and time. Press **OK** to confirm selection.

### TV connection

Select **TV Out** from the main menu and choose one of the existing settings. Press **OK** to confirm selection.

### Date and time

Select **Date/Time** from the main menu and set the current date and time. Press **OK** to confirm entry. You can also choose one of the following date formats: YY/MM/DD, DD/MM/YY or MM/DD/YY.

### Format SD card

Select **Format SD** from the main menu to format the installed memory card.

### Settings reset

Select **Sys. Reset** from the main menu to revert back to factory settings.

### Hardware version

Select **Version** from the main menu to check the current hardware version.

### Connecting the microscope to a TV-set

Plug the RCA cable into the microscope and connect it to your TV-set (refer to the User Guide of your TV-set for additional information). Turn on the TV and select the correct channel from the corresponding menu of the TV-set (refer to the User Guide of your TV-set for additional information).

## PortableCapture software

Insert the installation CD into your CD drive. The Main Menu window will be displayed. Select **PortableCapture software** to install the PortableCapture application. Keep in mind that a system restart will be required during installation. The installation CD contains a copy of this User Guide, which you can read by clicking the corresponding entry in the Main Menu.

Connect the microscope to your PC with a USB cable and run the PortableCapture application. A pop-up message will warn you if the microscope is not connected to the PC.

### Image resolution

From the **Options** menu, select **Preview Size** to set the required image resolution.

### Rotating the image

From the **Options** menu, select **Rotate** and choose to rotate created images 90° clockwise or counterclockwise.

### Date and time

From the **Options** menu, select **Date/Time** to set the date and time during preview.

### Language

From the **Options** menu, select **Language** to set the language.

### Full screen mode

There are three ways to enter full screen mode:

1. Double-click an image in the main window;
2. Click the **Full screen mode** icon on the toolbar;
3. From the **Options** menu, select **Fullscreen**.

To exit full screen mode, click the corresponding icon on the toolbar.

### Capturing an image

Click the corresponding icon on the toolbar or press the **F11** key on your keyboard. A thumbnail of the captured image will appear in the main window.

### Recording a video

Click the corresponding icon on the toolbar or press the **F12** key on your keyboard. The camcorder icon on the toolbar will change to the video recording icon. You can either click the corresponding icon on the toolbar or from the **Capture** menu.

*Note: recommended video resolution is 640x480.*

### Preview

Left-click the thumbnail of the image you wish to preview in the application. The image resolution will be displayed next to the thumbnail. Double-clicking a thumbnail will open the image in the main window.

### Saving images

To save an image, right-click the thumbnail of the image you wish to save. A drop-down menu will appear. Images may be saved in .jpg or .jpeg format, clicking **Advanced** allows you to save in other formats.

### Deleting images

To delete an image, right-click the thumbnail of the image you wish to delete. A drop-down menu will appear. When viewing images in a sequence, click the **Delete** icon on the toolbar to delete the image.

**Copying images**

To copy an image, right-click the thumbnail of the desired image and select **Copy** from the drop-down menu.

**Working with video clips**

Right-click a video clip thumbnail and choose whether you want to watch, copy or delete the chosen clip from a drop-down menu.

**Exiting the application**

To exit the application, choose **Exit** from the **File** menu.

**Calibration and measurements**

Place a ruler on the stage, focus your view and aimage. Take note of the current magnification by looking at the calibration scale on the microscope body. Double-click the image thumbnail to open it in a separate window. Enter the magnification value into the corresponding field in the upper right corner of the preview window. Upon calibrating the microscope, you will be able to measure various parameters of observed specimens. All the buttons on the toolbar of the preview window are described below.

**Line**

Press and hold the left mouse button to start drawing a line. Release the left mouse button to complete the line.

**Free-form line**

Press and hold the left mouse button and draw a line. Release the left mouse button to complete the segment. You can continue adding more segments to the free-form line.

**Radius**

Draw a straight line on the image. The application will automatically calculate the radius, as well as circumference and area of the corresponding circle.

**Diameter**

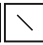
Draw a straight line on the image. The application will automatically calculate the diameter, as well as circumference and area of the corresponding circle.

**Angle**

Press and hold the left mouse button to draw a line on the image. Release the left mouse button and draw another line. The application will automatically calculate value of the resulting angle.

**Measurement units**

From the drop-down menu on the toolbar you can select measurement units for calculations (pixels, inches, kilometers, meters, centimeters, millimeters, microns).

Press this button to create

Press this button to edit fo

These buttons allow you to

Press this button to choos

Press this button to choos

Press this button to choos

Press this button to undo t

*Note:*  
1. *Magnification range of Levenhuk DTX : magnification values that exceed this range*  
2. *The application calculates values based on the magnification value entered. The results may not be precise.*  
3. *Measurement functions are only available for images.*

## Working with the application for Mac OS users

### Installing software

Browse to the **mac** folder on the installation CD and copy **PortableCapture** file to the desktop.

### Operating the microscope

Use a USB cable to connect the microscope to your computer. Run the PortableCapture application by double-clicking its icon.

### Observing specimens

A magnified image of the observed specimen may be seen in the main window of the application.

### Image resolution

From the **Options** menu, select **Preview Size** to set the image resolution. Keep in mind that the preview window cannot be resized (standard resolution is 640x480). The image resolution is shown in the bottom left corner of the preview window.

### Date and time

From the **Options** menu, select **Date/Time** to display or hide the date and time of capture during preview.

### Capturing an image

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu to capture an image. An image thumbnail will appear on the right side of the main window.

### Recording a video

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu to start recording a video.

*Note: during recording, the camcorder icon on the toolbar will turn red. Click the icon again to stop recording.*

### Editing images and videos

Image and video thumbnails are displayed on the right side of the main window of the application. Double-click an image thumbnail to open it in a separate preview window. Editing images with PortableCapture application is similar to working with other applications in Mac OS.

## Care and maintenance

Never, under any circumstances, look directly at the laser through this device, as this may cause PERMANENT EYE DAMAGE or BLINDNESS. White LEDs are very bright; do not look directly at them. Do not damage to your vision. Take necessary precautions when using the device. Do not use the device if you are wearing contact lenses or others who have not read or who do not fully understand the instructions. Do not use your microscope and before using it for the first time, read the instructions for every component and connection. Do not try to repair or modify the device for any reason. For repairs and cleaning of any kind, please contact the manufacturer. Do not center. Protect the device from sudden impact or vibration. Do not use excessive pressure when adjusting focus. Do not touch or clean the optical surfaces with your fingers. To clean the device, use only soft, lint-free wipes and special optics cleaning tools from Leica. Do not use the device for lengthy periods of time, and do not leave the device away from water and high humidity. Do not use the device to replace the dust cover after you are finished with the device. Do not use the device and stains. If you are not using your microscope, store the objective lenses and eyepieces separately from the device. Do not place away from hazardous acids and other chemicals. Do not use the device in other sources of high temperatures. To avoid damage to the device, turn off the application first, and then disconnect the microscope from the power illumination goes off completely, and disconnect the device from the power immediately if a small part or a battery is swallowed.

## Battery safety instructions

Always purchase the correct size and grade of batteries for the device. Always replace the whole set of batteries at one time. Do not mix old ones, or batteries of different types. Clean the device before and after battery installation. Make sure the battery polarity (+ and -). Remove batteries from equipment after use. Do not use the device for a long period of time. Remove used batteries promptly after use, as this may cause leakage, fire, or explosion. Do not use the device in high temperatures, leakage, or explosion. Never use the device. Remember to switch off devices after use. Keep the device away from the risk of ingestion, suffocation, or poisoning.

## Specifications

Image sensor	5Mpx (up to 12Mpx by interpolation)
Magnification	20x–200x–500x
Available image resolutions	12Mpx, 8Mpx, 5Mpx, 3Mpx, 1.3Mpx, VGA
Photo/video format	*.jpeg/*.avi
Memory card	microSD up to 32Gb (not included)
Objective lens	high-quality lens
Display	color display 3" TFT, 4:3 ratio
Illumination	8-LED system with variable brightness
Power source	Li-ion battery (3.7V/1050mAh)
	Operation time: 4 hours, charging time: 5 hours
Frame rate	30fps
Power cable	100–240V, 50/60Hz; 5V, 1A output
OS language	English, German, French, Spanish, Italian, Portuguese, Japanese, Chinese
Software language	English, German, French, Spanish, Russian, Italian, Portuguese, Polish, Dutch, Japanese, Korean, Chinese
Dimensions (HxWxD)	330x70x150mm
Weight	700g (with the battery)

Levenhuk reserves the right to modify or discontinue any product without prior notice.

## System requirements

Operating system: Windows XP SP2/Vista/7/8, Mac 10.6–10.8,  
CPU: at least P4 1.8GHz, RAM: 512MB, interface: USB 2.0

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a lifetime warranty against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk. This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

For further details, please visit our web site: [www.levenhuk.com/warranty](http://www.levenhuk.com/warranty)

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Levenhuk® is a registered trademark of Levenhuk, Inc. © 2006-2015 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 924-D East 124th Ave. Tampa, FL 33612 USA