# Levenhuk biological microscopes

Levenhuk 850B (binocular)

Levenhuk 870T (trinocular)

Levenhuk D870T (trinocular)

User Manual

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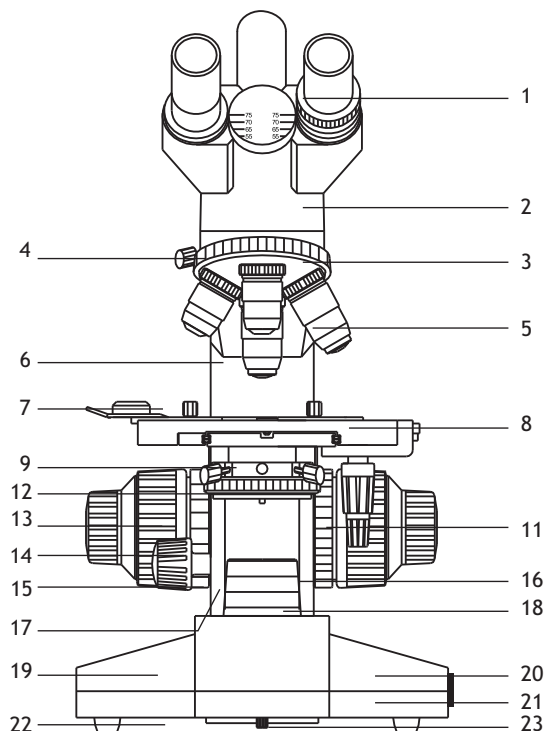
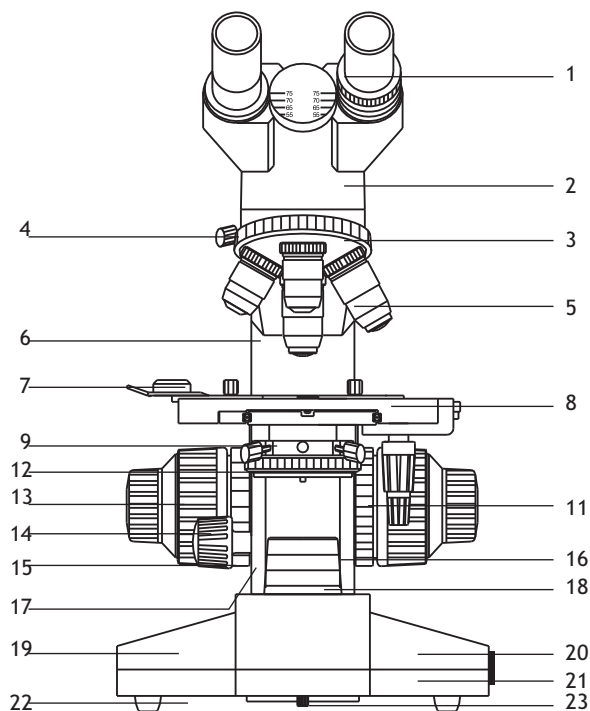
**levenhuk**<sup>®</sup>  
Zoom&Joy

- 1 Eyepiece
- 2 Binocular head / Trinocular head
- 3 Revolving nosepiece
- 4 Locking screw
- 5 Objective lens
- 6 Arm
- 7 Specimen holder
- 8 Stage
- 9 Abbe condenser
- 10 Stage left/right movement control
- 11 Iris diaphragm
- 12 Fine focus knob
- 13 Coarse focus knob
- 14 Condenser vertical movement control
- 15 Focus system tension adjustment ring
- 16 Stage locking knob
- 17 Collector lens
- 18 Base
- 19 Power switch
- 20 Brightness adjustment knob
- 21 Rubber feet
- 22 Lamp compartment locking screw

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**Never look directly at the Sun through optical devices or with the unaided eye! It can result in permanent eye damage or even BLINDNESS.**

## General Information

These microscopes are high precision instruments that may be used for clinical research and university studies; in laboratories and hospitals, technological centers and research institutes. They also perform well in medical studies in biology, bacteriology, cytology and pharmacology. Levenhuk 850B and 870T may be used for microphotography with additional accessories. The kit of Levenhuk D870T already includes an eyepiece digital camera that allows you to observe specimens on your PC monitor and save images on the hard drive.

This manual contains detailed information on the design and operational principles of your new microscope, as well as a troubleshooting and care and maintenance section. It is highly recommended that you read this manual before using the microscope for the first time. Abiding by the following instructions will ensure maximum service life of your microscope:

- This instrument was designed for high-power observations of various biological specimens. It should not be used for any other purpose.
- Do not try to take this microscope apart on your own, as you might damage the intricate optical system within. If the troubleshooting section of this manual did not contain a solution to your issue, you should contact the local Levenhuk service center.
- Always unplug the microscope from a power source before opening the base or changing the illumination lamp. Regardless of the lamp type (halogen or incandescent), give it some time to cool down before trying to change it, and always change it to a lamp of the same type. When using the microscope, try not to use it near flammable materials or substances (benzene, paper, cardboard, plastic, etc.), as the base may heat up during use, and might become a fire hazard.
- Always use the power supply with the proper voltage, i.e. indicated in the specifications of your new microscope. Plugging the instrument into a different power outlet may damage the electric circuitry of the microscope, burn out the lamp, or even cause a short circuit.
- Never touch the microscope optics with your fingers, as fingerprints and smudges will significantly reduce the quality of your views.
- These microscopes perform best under the following conditions: room temperature from 0 °C to 40 °C (32-104 °F), relative humidity not exceeding 85%. High temperatures and humidity may cause lenses to fog from inside, thus damaging the optical system.
- Any microscope is a very delicate instrument, so it should be handled with care. Even a slight impact might damage it.

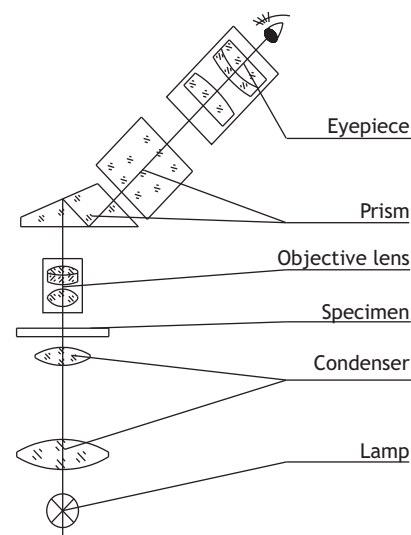
## Operating principle

Light emitted from a lamp passes through the condenser and illuminates a specimen. The image is magnified by the objective lens, refracted by the prism and then magnified once more by the eyepiece.

Total magnification = (objective lens magnification) x (eyepiece magnification).

The kit of these microscopes includes two WF10x eyepieces with wide field of view:

Magnification	10x
Field of view	18 mm (0.71 in)
Working distance	24.95 mm (0.98 in)

## Objective lenses

The standard kit of these microscopes includes the following DIN achromatic objective lenses: 4x, 10x, 40x (spring-loaded) and 100x (spring-loaded, oil immersion). When using the 100x objective lens, both the lens and the cover glass of the observed specimen should be immersed in a transparent oil of high refractive index. This displaces the air and, as a result, increases the quality of your views. If some air bubbles are still present, you may have to rotate the revolving nosepiece once, or add some more immersion oil. Upon finishing the high-power observations, you should wipe the objective lens and the cover glass clean. Otherwise, the oil will dry and reduce the quality of subsequent views.

DIN achromatic objective lens	4X	10X	40X	100X
Magnification	4	10	40	100
Aperture	0.1 mm	0.25 mm	0.65 mm	1.25 mm
Working distance	36.9 mm (1.45 in)	7.116 mm (0.28 in)	0.632 mm (0.025 in)	0.158 mm (0.006 in)
Comments			Spring-loaded	Spring-loaded, oil immersion

## Binocular head

The binocular head of these microscopes may be inclined at a 30° angle and rotated 360 degrees. The interpupillary distance adjustment range is 55-75 mm (2.16-2.95 in), the diopter adjustment range is from +5 to -5 diopters on each eyepiece.

## Revolving nosepiece

The state of the art nosepiece provides for smooth rotation and easily locks in place on the desired objective lens. The parfocal objective lenses used in these microscopes provide for simple switching from one magnification to the other, without having to refocus your view over and over again.

## Stage

These microscopes are equipped with a 140 mm x 155 mm (5.5 in x 6.1 in) stage with coaxial controls. Press the stage clip lever and place a slide with a specimen on the stage. Release the lever and it will securely lock the slide in place. You can move the stage horizontally by using the coaxial controls below the stage. It moves within 70 mm (2.76 in) range in 0.1 mm (0.004 in) increments.

## Abbe condenser

The numerical aperture (N.A.) of this high quality condenser is 1.25. Its position may be adjusted by turning the locking screws, without any additional tools:

- rotate the revolving nosepiece until either a 4x or a 10x objective lens is locked in place;
- rotate the iris diaphragm to reduce its diameter;
- lower the condenser until the diaphragm image is somewhat sharp;
- focus the view by loosening or tightening the locking screws.

The condenser has been aligned with the objective lens during production. If this alignment hasn't been performed, the field of view will be very irregular, with parts of it very bright, and others very dark. The rack-and-pinion mechanism moves the condenser up and down within a 20 mm (0.79 in) range. Generally, you would want to raise the condenser when using high-power objective lenses (40x, 100x) and lower it with low-power objective lenses (4x, 10x). The highest position of the condenser is locked at 0.2 mm away from the stage.

If the condenser-objective lens alignment has somehow gone awry, you have to perform the following operations: loosen the nut on the locking screw;

- place a slide with a specimen on the stage;
- raise the condenser to its highest position;
- tighten the locking screw until the condenser touches the bottom of the stage;
- retighten the nut on the locking screw.

Rotate the iris diaphragm to adjust its diameter in 2-30 mm (0.08-1.18 in) range, so that it corresponds with the N.A. of the objective lens currently in use. When the iris diameter is approximately 70-80% of the objective lens N.A., the observed image is sharp and clear. Remove one of the eyepieces and look through the tube. You will see the iris diaphragm. You can use a color filter with the diaphragm for additional contrast and detail on the image. The filters come in blue, green or yellow.

## Focus system

The coarse and fine focus systems are operated with a rack-and-pinion mechanism. The coaxial design of the focus system allows you to adjust coarse and fine focus without any difficulties. Fine focus system adjusts in 0.002 mm (0.00008 in) increments; total adjustment range is 25 mm (0.98 in). You can use the coarse focus knob to quickly raise or lower the stage, and fine-tune the view with the fine focus knob.

Upon moving the stage into a desired position, you can secure it in place with the stage locking screw. Rotate it clockwise to lock the stage in its current position. It will remain in this position until you loosen the locking screw. If you are observing a large and heavy specimen, you might want to increase the coarse focus system knob tension. Rotate the adjustment ring on the focus knob and the stage will be lowered slowly and steadily, even with heavy specimens.

## Illumination

Generally, a 6V/20W halogen lamp is used in the illumination system of a microscope. To adjust its brightness, rotate the corresponding knob, to better illuminate the observed specimen. The illumination should not be too bright with low-powered objective lenses and vice versa. When changing the lamp to a new one, make sure to align it along the optical axis, otherwise, the quality of your views will be significantly decreased.

To change the old lamp, loosen the screw on the bottom of the base, remove the cover and replace the lamp.

## Assembly

Unpack your new microscope with great care. Check if all the parts and accessories are present. Note that some parts might have been disassembled for easier shipping, therefore, you will have to assemble them before you can begin your observations.

- Assemble the binocular head. Loosen the locking screw and remove the plastic cover from the arm. Remove the plastic cover from the binocular head. Attach the binocular head to the arm and secure it in place with a locking screw. Do not try to loosen the other two screws on the arm.
- Assemble objective lenses. Usually, we assemble the objective lenses on the factory lines, but in some models, they are shipped separately. First, remove the dust cap from the nosepiece and take out the objective lenses from their packaging. Attach the objective lenses to the nosepiece in ascending order (from lowest magnification to highest).
- Assemble eyepieces. Remove the dust cap from the tube and attach the eyepieces.

## Using the microscope

- Choose a spot, where the microscope will not be in direct sunlight. Try to move as far away from windows as possible and do not point eyepieces toward windows. Direct sunlight may damage the optics and lower the contrast of images.
- Place a slide with a specimen on the stage and lock it in place. Make sure that the cover glass faces upwards, otherwise, you might have problems focusing your view at higher magnifications (40x or 100x).
- Use a low-power objective lens (4x or 10x) and turn the illumination on.
- Using the coaxial stage movement controls, position the slide, so that the specimen is centered above the condenser lens.
- Focus your view with the coarse focus knob. It is much easier to focus your view at lower magnifications, so that when you use a high-power objective lens, you will only have to fine-tune the focus with the fine focus knob.
- Now that your view is focused, rotate the revolving nosepiece to a high-power objective lens. Thanks to the parfocal lenses in these microscopes, you will only have to use the fine focus knob to bring your views into focus. It has to be mentioned that 4x and 10x objective lenses will never touch the surface of the slide, due to the built-in limiters. And although 40x and 100x may reach the slide, thanks to their spring-loaded construction, they will never break it.
- If you cannot focus your view, you might want to try changing the diameter of the iris diaphragm, so that it corresponds to the N.A. of the objective lens used.

## Troubleshooting

If you are experiencing some problems with your instrument during its service, refer to this section for possible solutions. However, if this doesn't help, you should contact your local service center. Remember, do not try to take the microscope apart, as this will make your warranty null and void, as well as may lead to irreversible damage to the intricate optical system of the instrument.

### Problems with the microscope:

Problem	Possible cause	Solution
Impossible to focus the view	The stage is too low	Raise the stage with the coarse focus knob
The objective lens often damages slides	The stage is too high	Lower the stage with the fine focus knob
Impossible to focus the view when using high-power objective lenses	The slide may have been placed upside down, or the cover glass is too thick	Place the slide properly or opt for a standard cover glass (0.17 mm (0.007 in) thick)
When rotating the nosepiece, objective lenses always scrape the surface	The cover glass is too thick	Opt for a standard cover glass (0.17 mm (0.007 in) thick)
Impossible to move the slide smoothly	The stage clips are not locked	Lock the stage clips manually
Incomplete binocular vision	Incorrect interpupillary distance	Adjust the interpupillary distance
	Incomplete diopter adjustment	Readjust diopter settings
	Insufficient brightness	Check the illumination source and readjust the brightness

### Problems with the optical system:

Problem	Possible cause	Solution
The field of view is reduced or not fully illuminated	The objective lens is not locked in place	Rotate the revolving nosepiece to lock the desired lens in place
	The optical axis is not aligned	Readjust the illumination lamp
	The optics of the objective lens is dirty or dusty	Clean the objective lens
The field of view is reduced or not fully illuminated	The optics of the objective lens is dusty or dirty	Clean the objective lens
	The surface of the slide is dusty or dirty	Clean the slide
	The condenser lens is too low	Raise the condenser lens
Poor quality of views: reduced contrast, blurred details, etc.	The specimen is not protected with cover glass	Put the cover glass over the specimen
	The cover glass is too thick or too thin	Opt for a standard cover glass (0.17 mm (0.007 in) thick)
	The slide is placed upside down	Place the slide correctly
	The upper lens of the objective lens is dirty	Clean the lens
	The optics of the objective lens is dirty	Clean the objective lens
	Immersion oil was not used at high magnifications	Add immersion oil
	Air bubbles are present in the immersion oil	Rotate the revolving nosepiece or add some more oil

Poor quality of views: reduced contrast, blurred details, etc.	Using immersion oil with low refractive index	Use the required immersion oil with high refractive indices
	Prism surface is dirty	Clean the prism
	Iris diaphragm diameter is too wide or too small	Adjust the diameter
	Condenser lens is too low	Raise the condenser lens
One side of the field of view is darker than the other	Condenser lens is placed incorrectly or is inclined.	Adjust the condenser lens
	The objective lens is not aligned to the optical axis	Rotate the revolving nosepiece until the objective lens locks in place
The view is constantly shifting during observations	The slide is not locked on the stage	The objective lens is not aligned to the optical axis
	The objective lens is not aligned to the optical axis	Rotate the revolving nosepiece until the objective lens locks in place
The observed image is yellow	A blue filter is not used	Use the blue filter
Condenser lens is too low	Iris diaphragm diameter is too low	Adjust the diameter
	Condenser lens is too low	Raise the condenser lens
	The optics of the objective lens is dirty	Clean the objective lens

#### Problems with electric parts:

Problem	Possible cause	Solution
Illumination does not work, even though it is turned on	Power cable has disconnected	Check the power cable
	The lamp is not installed	Install the lamp
	The fuse is blown	Replace the lamp or the fuse
Lamps burn out quickly	Incorrect lamp is being used	Use the standard lamp
	Lamp voltage is too high	Lamp voltage is too high
The illumination is too dim	Incorrect lamp is being used	Use the standard lamp
	The power source voltage is too low	Lamp voltage is too high
The lamp keeps blinking, brightness keeps changing	Electrical surges	Adjust the power source
	The filament is about to burn out	Change the lamp
	Power cable has disconnected	Check the power cable

## Care and maintenance

- Unpack the microscope with great care, make sure nothing falls on the ground;
- All the optics are calibrated on the factory line, do not try to take them apart on your own;
- The binocular head and the focus system are high-precision elements, and should not be taken apart. If you have any problems with either of them, contact our local service center;
- All mechanical elements have to be protected from dust; all the mechanisms should be greased regularly; all the optical elements should be kept clean;
- Store your microscope in a clean, dry place. Upon finishing your observations, replace the dust cover and unplug the instrument from the power source.

## Digital camera

Levenhuk D870T comes with an 8 Mpx C800 NG digital camera.

The camera allows you to observe specimens in fine detail and true colors on your PC monitor and save images on the hard drive.

The special software that comes in the kit allows you to view and edit the resulting images. Supported file formats include: \*.bmp, \*.jpg, \*.jpeg, \*.png, \*.tif, \*.tiff, \*.gif, \*.psd, \*.ico, \*.emf, \*.wmf, etc.

Connect the camera to your PC via a USB cable that also acts as a power cable, so you do not have to plug it into a power source.

Compatible OS: C800 NG - Windows XP/Vista/7/8.

Camera model	C800 NG
Max. Resolution (for still images)	3264x2448
Megapixels	8.0
Sensor	1/2.2" CMOS
Dimensions (cylind unit maximum diameter)	50 mm (1.97 ")
Pixel size	1.67µm x 1.67µm
Sensitivity, V/lux.sec@550 nm	0.31
Mounting location	Eyepiece tube (replaces the eyepiece)

Image format	BMP, TIFF, JPG, PICT, SFTL, etc.
Field of view diameter	18 mm (0.71 in)
Spectral range	400 nm - 650 nm
Exposure	ERS
White balance	Auto/manual
Exposure control	Auto/manual
Software image control	Image size, brightness, exposure control
Interface	USB 2.0, 480 Mb/s
System requirements	Windows XP/Vista/7/8; USB port
Software	Software for image viewing and editing
Shell	Metal cylinder, painted black
Power source	USB 2.0 cable
Temperature range	-30 - 70 °C / -22-158 °F

Generally, when using a low-power objective lens (4x, 10x) the image on the PC monitor will be too bright. In this case you can:

- reduce the illumination brightness;
- reduce the iris diaphragm diameter;
- lower the Abbe condenser lens.

When using a high-power objective lens (40x, 100x) the image on the PC monitor will be too dark. In this case you can:

- increase the illumination brightness;
- increase the iris diaphragm diameter;
- raise the Abbe condenser lens.

## Caution

**Never look at the sources of bright light or lasers through the camera or the microscope: it will cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS!**

**Do not try to take the camera or the microscope apart on your own.**

**Protect the camera and the microscope from moisture; do not use them in rainy weather.**

**Protect the camera and the microscope from impacts or excessive pressure.**

**Do not overtighten the locking screws.**

**Keep the camera and the microscope away from hazardous environments, home and care heaters, incandescent lamps or open fire.**

**When cleaning the lenses, first blow any dust or debris off the surface or wipe the off with a soft brush. Then wipe the lens with a soft cleaning cloth slightly damp with spirits or ether.**

**Never touch the optics with your fingers.**

## International Warranty

- The Vendor guarantees that the quality of the Levenhuk product purchased complies with the technical documentation requirements on conditions that the consumer observes rules of transportation, storage conditions and operating instructions.
- Levenhuk Ltd. warrants the product against defects in materials.
- During the warranty period the Buyer can return the defect product to the Vendor or to Levenhuk service center. Levenhuk Ltd. will repair or replace the defect product at its own discretion.
- No claims are accepted in case the properly filled warranty slip is absent or contains corrections, or if the defect product has not been provided by the Buyer. The manufacturer or the seller is not liable for any damages caused by misuse of the product. No guarantee is provided in cases the product was used for purposes other than that intended or has mechanical damages, scratches, cracks, optics damages, as well as in cases the product is broken-down due to impact damages, squeezing or stretching or the product has been disassembled or repaired by unauthorized personnel.
- This warranty does not cover components and/or accessories with limited use period, batteries, etc.

The period of warranty is 3 (three) years beginning on the date of purchase. Please keep the warranty slip along with your receipt.

For any warranty services please contact the local Levenhuk branch.

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Levenhuk® is a registered trademark of Levenhuk, Ltd.
Copyright © 2006-2013 Levenhuk Ltd. All rights reserved
Developed and manufactured for Levenhuk, Ltd.